

# Instruction manual

## *For safety cabinet*



### Model ASVM11

## CONTENT

---

|            |                                |        |
|------------|--------------------------------|--------|
| <b>1.</b>  | Description / Use              | Page 2 |
| <b>2.</b>  | Technical features             | Page 2 |
| <b>3.</b>  | Installation and commissioning | Page 2 |
| <b>4.</b>  | Internal equipments            | Page 3 |
| <b>5.</b>  | Storage                        | Page 3 |
| <b>6.</b>  | Ventilation                    | Page 3 |
| <b>7.</b>  | Safety controls                | Page 4 |
| <b>8.</b>  | Spare parts                    | Page 5 |
| <b>9.</b>  | Recycling                      | Page 5 |
| <b>10.</b> | Warranty                       | Page 5 |
| <b>11.</b> | Control sheet for the user     | Page 7 |

## 1. DESCRIPTION / USE

The chemicals have to be stored in safety cabinets. These safety cabinets allow storing either flammable products, acids, bases or toxics (low dangerous), or several types together.

## 2. TECHNICAL FEATURES

### Compliance :

- Compliance with INRS recommendations
- Compliance with the articles R. 5132-66 and R. 5132-68 of the Public Health Code
- Standardized pictograms compliant with the standards ISO 3864, ISO 7010, NF X8.003 and the directive 92/58/CEE

### Technical features :

- Construction in 10/10th steel
- White epoxy paint RAL 9010
- Standardized pictograms provided 1x « flammables », 2x « corrosives » and 1x « toxics »
- Manual doors
- Locking system
- Shelves adjustable in height, every 100 mm
- Air inlet and outlet
- External dimensions : H 1825 x W 1100 x D 545 mm
- Storage capacity : 260 liters (2x 130)
- Retention capacity : 92 liters (2x 46L)
- Loading capacity : 50 KG per shelf
- Weight : 115 KG

## 3. INSTALLATION AND COMMISSIONING

### Installation

The safety cabinet has to be installed on a flat surface that can bear its weight with the products inside.

Ensure that the safety cabinet is correctly wedged, as for the cabinet not to topple. The eventual blocking has to be done by the user.

The correct leveling of the cabinet must cart off any distortion of the frame, allow the doors to be strictly parallel compared to the frame, and take away any rubbing of the doors on the cabinet's frame.

The installation location must guarantee :

- That it is not exposed to any heat source
- That it is protected against moisture and that the ambient temperature of the location is neither less than 5°C, nor more than 40°C

### Transport and handling

The safety cabinet that you received has been protected and transported in optimal conditions, as to be protected from any damage.

It has to be lifted only by suitable equipment (suction cup, strap...). It has to be carried in a vertical position, and it must not be lying. It is imperative not to tilt the cabinet when it is carried. Indeed it can cause damages that can alter the right functioning of the cabinet.

On the other hand, if it has to be moved, the cabinet has to be placed on a pallet and strapped, as to avoid any fall.

### **Commissioning**

After having unwrapped the cabinet, ensure that you also received a set of two keys.

Your safety cabinet has to be installed, used and maintained as to guarantee the users protection.

## **4. INTERNAL EQUIPMENTS**

The cabinet is provided with 6 retention shelves and 2 retention tanks. The retention capacity per shelf is 7 liters and per retention tank 25 liters.

Take care not to exceed the loading capacity per shelf of 50 KG.

The shelves must be fixed at the chosen height thanks to the rack and pinion.

## **5. STORAGE**

Your safety cabinet is exclusively destined to store low risk chemicals in laboratories, industries, schools... in closed containers.

It must not store high concentrated corrosive products or very flammable products.

We remind you that the storage of corrosives can alter the quality of your safety equipment. On the other hand, we remind you that damages due to a bad use of the equipment and in particular to corrosion won't be covered by our warranty.

## **6. VENTILATION**


We highly recommend connecting your safety cabinet to a ventilation system, as to evacuate the eventual vapors contained in it.

The air renewal has to be at least 10 times the volume of the cabinet, per hour. Moreover, ensure that the ventilation runs 24h/7.

### **Connection to a ventilation system**

The air inlets are situated at the bottom, on the doors ; the outlet is situated on the top. See the schema below.

|                                                        |                                                                                                                                                                                                                                |
|--------------------------------------------------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| <b>Model</b><br><b>ASVM11</b>                          | <b>ECOSAFE</b><br>www.ecosafesa.com                                                                                                                                                                                            |
| Scale 1 : 15<br>Measure unit : mm<br>Date : 06/03/2019 | Chemin des Champs Courbes 15 –<br>Zone Industrielle de Vallaire Est<br>CH-5204 Ecublens<br>Tél : +41 (0)21 694 10 04<br>ECOSAFE reserves the right to<br>discontinue or change specifications of<br>any device without notice. |

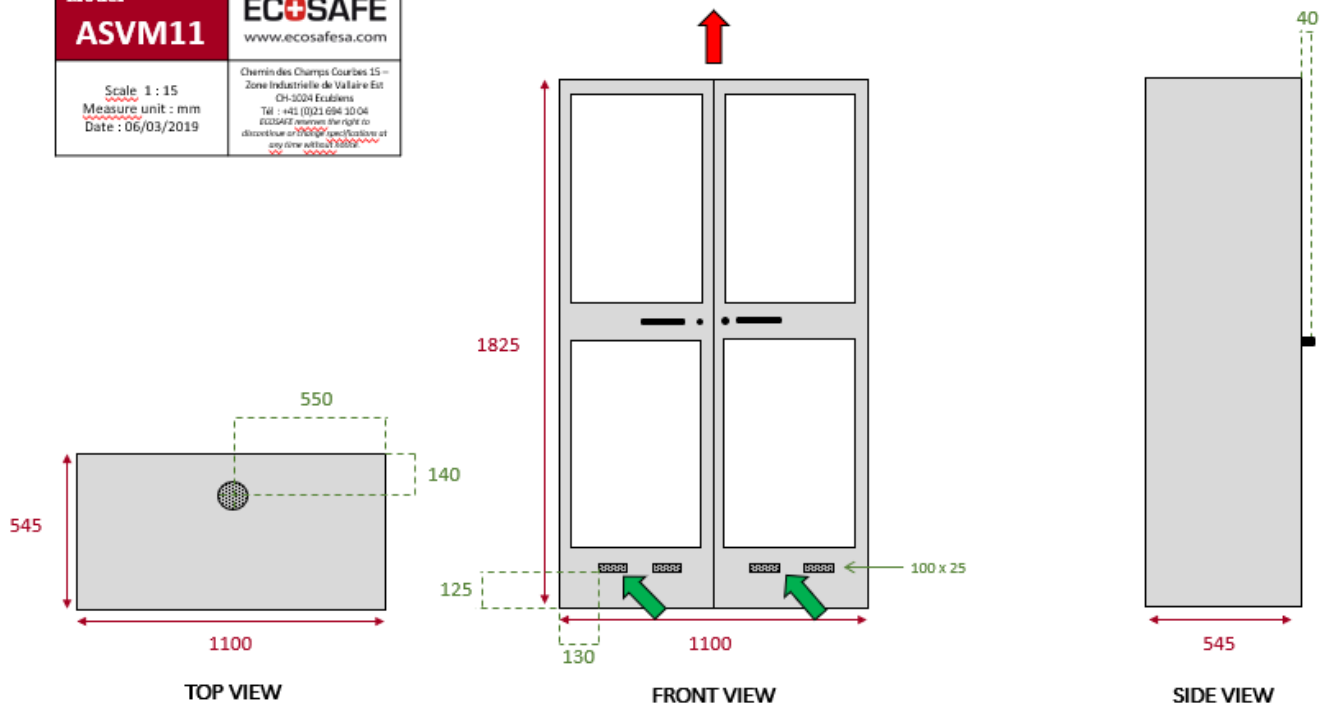

### How to choose a ventilation or a filtration box ?

**Ventilation boxes** are connected to the safety cabinet with a connecting kit (including a connecting sleeve, 2 hose clamps and 1 meter of flexible sheath) ; the fan included in the box extracts the toxic vapors contained in the cabinet, and reject it to the outside of the building thanks to the roof or wall extraction.

If you cannot reject the vapors to the outside, you need a filtration box.

**Filtration boxes** are equipped with an activated carbon filter, which purifies the toxic vapors in the box, and rejects it in the room in a healthy fashion. Filtration boxes are delivered with the activated carbon filter that you need to renew once a year.

### Running without any mechanic ventilation

It is possible for your safety cabinet to run without mechanical ventilation. However, your cabinet has to be placed in a location that provides its own ventilation system (at least 5 times the volume of the location, per hour).

The cabinet itself and a radius of 2,5 meters directly around is a dangerous area subject to explosions.

On the other hand, if the ventilation system of your cabinet is not running, you have to indicate it, by putting an etiquette on the cabinet, saying "Safety cabinet not connected to the mechanic ventilation" (etiquette is not provided).

## 7. SAFETY CONTROLS

Your safety cabinet is a safety equipment. As to guarantee its perfect efficiency, you have to check the equipment regularly.

**Daily safety controls :**

- Check if the products (liquids, powders...) have flowed in the containment sumps. If it has, they have to be emptied and cleaned promptly, with suitable equipment (absorbent mat...).
- Check that the cabinet and its internal equipments are not corroded.
- Check that the cabinet does not have any mechanical damages.
- Check the proper closing of the doors.
- Check the proper working of the locking system.

**Monthly safety controls :**

- Check the proper working of the doors (opening and closing)
- Check the working of your ventilation system thanks to, for example, a piece of paper  
(*Only if your cabinet is connected to a mechanical ventilation system*)

**8. SPARE PARTS****Elements :**

- Handle : *Contact us*
- Keys : *Contact us*
- Standardized pictogram : P/N **PICTO**
- Additional shelf : P/N **EE2**
- Protective shelf in PVC : P/N **SEPVC**
- Retention tank in PVC : P/N **SBPVC**

**9. RECYCLING**

The safety cabinet can be entirely dismantled. Various elements such as the steel can be recycled. Ensure to respect the national and local waste disposal terms. Those elements should not be mixed with household waste.

**10. WARRANTY**

Your equipment holds a 12 months warranty against every manufacturing defect (limited to ordinary conditions of use, from the date indicated on the delivery note).

Our warranty can be applied only in case of manufacturing defects or hidden fault. The warranty covers only the replacement of the faulty goods or pieces indicated by our quality or technical department, as they are the only one able to determine the condition. Repair of the equipment must be performed in our warehouse.

The spare parts are under 3 months warranty upon their settlement.

The warranty cannot be applied in the following cases :

The equipment has not been stored, used or maintained in accordance with the standards by the user

In case of a damage of the equipment by the user or a third party

The equipment had a modification or a repair by the user or a third party, without any prior written authorization by ECOSAFE

In case of a bad use of the equipment

In case of a normal attrition of the equipment

In case of a bad installation of the equipment

In case of damage caused by the corrosion

The activation of the warranty won't extend the term. Our warranty stops if the customer has not notified us in a 7 days term after seeing the fault ; date that the customer has to prove.

